

PHARMA MACHINERIES

COMMUNITING MILL



MINI COMMUNITING MILL





RIDDHI PHARMA MACHINERY LTD.

CIN: U33112GJ1995PLC028504

Manufacturer & Exporters of Pharmaceutical Machineries

MUMBAI OFFICE: Om-shivam soc., Tarun Bharat Bldg. No.2, Opp. Cigaratte Factory, Sahar Road, Andheri (East), Mumbai - 400 099. Maharashtra, India Tel.: +91-22-28321332 / 28382479 Mobile: +91- 9821018579 / 9969994187 Email: <a href="mailto:hkshah@riddhipharma.com/riddhipharma@yahoo.com/riddhipharma Web: www.riddhipharma.com / www.riddhipharma.net

REGD. OFFICE & FACTORY: Riddhi Estate, Gulab Nagar, Opp. AEC, Amraiwadi, Ahmedabad - 380026. India.





TYPE OF SCREENS AVAILABLE:

TYPE 'A' SCREENS - Perfor

Perforated Round Holes: 3/32", 7/64", 1/8", 5/32", 3/16", 7/32", 1/4", 5/16", 3/8", 1/2", 5/8", 3/4" and 1". Perforated Square Holes: 1/4", 3/8", 1/2", 7/8" and 1 1/8".

TYPE 'B' SCREENS -

Perforated Round Holes with back reinforcing plate: 1/50", 1/42", 1/37", 1/32", 1/25", 1/20" and 1/16".

TYPE 'C' SCREENS -

Woven Stainless Steel Wire Mesh with back reinforcing plate: 3, 6, 8, 10, 12, 14, 16, 20, 30, 40 and 50 mesh.

TYPE 'D' SCREENS -


Woven Stainless Steel Wire Mesh mounted on front and back reinforcing plate of 36 Square Holes: 60, 80 mesh.

TYPE 'E' SCREENS -

Woven Stainless Steel Wire Mesh mounted on front and back reinforcing plate of 36 Square Holes: 100, 200Mesh

Perforated screens are also available with Holes size in mm.



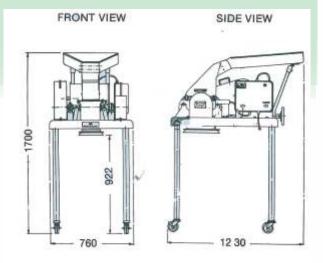

ALL DIMENSIONS ARE IN MM



RIDDHI PHARMA MACHINERY LTD.

Manufacturer & Exporters of Pharmaceutical Machineries

OFFICE: MUMBAI Om-Shivam Bldg.,2, Tarun Bharat, Sahar Rd, Andheri(E), Mumbai-400 099.Maharashtra, India. Tel.: 91-22-2832 1332 / 2838 9432 Fax: 91-22-2838 2479 Res: 91-22-2683 3777, Mob: 98210 18579

Whatsapp: 919821018579 Skype: hiten.k.shah

E-mail: hitenkshah@hotmail.com/ hitenks@vsnl.com **Website:** www.riddhipharma.com **FACTORY:** Riddhi Estate, Gulab Nagar, Opp. AEC, Amraiwadi, Ahemdabad-380 026. India.



RIDDHI MILL COMMINUTING MACHINE PLAIN & WATER JACKETTED MODEL: RDFM

<u>APPLICATION</u>: Riddhi mill has been specifically designed for wet and drygranulation, pulverization or dispersion of products oringredients found in the pharmaceautical, Chemical and foodindustries. It is also finds application in Cosmetics, confectionery, Fertilizers, Insecticides, detergents and Electroflux Industries etc.

<u>DESCRIPTION</u>:Latest improvements in design makes this unit a truely GMP model and Versatile in operation & maintenance.

Unit consist of S.S.304 quality hopper, feed throat and processing chamber (can be provided in S.S.316 quality at extra cost) with beaters assembly , heavy duty motor, star delta starter with overload relay, three speed step pulley, suitable screen , S.S. cladded base having tubular legs with castors,

optional jacket for the feed throat and processing chamber forcooling. <u>OPERATION</u>: Product kept in the hopper slide down to the enclosedchamber, where it is milled in between beaters and screento the required mesh/micron size which can be collected the bottom of the container. Quality and output depends upon the three main variables 1) Beaters shape (knife/impact forward) 2) Speed and

3)Screen size.Knife edge for coarse or large particles and impact or flatedge for finer particles. Highspeed for fine griding, lowspeed for coarser particles. A wide range of screens is available for various process needs.

TECHNICAL SPECIFICATION

OUTPUT : 50 kgs. to 800 kgs per hour depending on the product and the reduction required in size.

ROTOR SPEED : 1000/2250/4500 RPM.

BEATERS : 16 NOS Stainless steel with knife and impact edges

SCREEN AREA : 495 mm x 135 mm (66825sq.mm)

MOTOR : 7.5HP/1400RPM/440V/50CS/3PH/AC.

STARTER: Star Delta with no-load and overload protection.

NET WEIGHT : 400 kgs. GROSS WEIGHT : 530 kgS.

DIMENSIONS : 1230 x 760 x1575 MM(HT)49 x30 x62 inch app.

CASE DIMENSIONS: 1380 x1050 x1900 MM(HT)

55 x 42 x 75 inch



RIDDHI PHARMA MACHINERY LTD.

CIN: U33112GJ1995PLC028504

Manufacturer & Exporters of Pharmaceutical Machineries

MUMBAI OFFICE: Om-shivam soc., Tarun Bharat Bldg. No.2, Opp. Cigaratte Factory, Sahar Road, Andheri (East), Mumbai - 400 099. Maharashtra, India Tel.: +91-22-28321332 / 28382479 Mobile: +91- 9821018579 / 9969994187 Email: hkshah@riddhipharma.com /riddhipharma@yahoo.com

Web: www.riddhipharma.com / www.riddhipharma.net

REGD. OFFICE & FACTORY: Riddhi Estate, Gulab Nagar, Opp. AEC, Amraiwadi, Ahmedabad - 380026. India.

